

Technical Note: Comparison between SEC35 and SEC70 columns for extracellular vesicle isolation

Introduction

Size exclusion chromatography (SEC) is a widely used technique for the separation of extracellular vesicles (EVs), including exosomes and microvesicles, from biological fluids or culture media. At Immunostep, we offer two types of SEC columns: **SEC35** and **SEC70**, adapted to different experimental needs depending on particle size, desired purity, and sample type.

Main differences between SEC35 and SEC70

Feature	SEC35	SEC70
Resin pore size	Small	Large
Separation range	~35–350 nm	~70–1000 nm
Target particles	Exosomes, small EVs	Microvesicles, medium and large EVs
EV recovery	High	Moderate
Purity (free of proteins/lipoproteins)	Moderate	High
Protein co-elution	May contain some albumin or HDL	Lower presence of contaminants
Ideal applications	Cases where recovery and purity are equally important	When purity is a priority (e.g. proteomic or functional analysis)

Recommendations for use

- **SEC35:**
Ideal for researchers working with exosomes or small EVs who need **maximum particle recovery**, even if purity is slightly lower.
Common applications: RNA studies, NTA, subsequent immunocapture.
- **SEC70:**
Recommended when **purity is key**, especially in samples with high protein load (such as plasma). Effective for separating EVs from lipoproteins and albumin.
Common applications: proteomic analysis, functional studies, diagnostic use.

Additional considerations

- Both columns are compatible with samples such as **plasma, serum, urine, and culture media**.
- It is recommended to use fractions between **F7 and F12** to obtain EVs with greater integrity and less contamination.


Format and compatibility

- The columns are available in a ready-to-use format.
- They can be combined with immunocapture kits (such as ExoStep™) for subsequent phenotypic characterisation by **flow cytometry, ELISA, or surface labelling**.

Conclusion

The choice between SEC35 and SEC70 will depend on the desired balance between **recovery and purity**, as well as the size of the EVs to be analysed. For personalised assistance, the Immunostep technical team is available to help you select the best option for your specific application.

Technical contact:

 tech@immunostep.com

 www.immunostep.com

Immunostep S.L. – Specialists in the detection and characterisation of extracellular vesicles